Instructions / Instrukties
Instruktionen / Instrucciones
Instruções

 **BRADY®**

www.bradylabid.com

While every precaution has been taken in the preparation of this document, Brady assumes no liability to any party for any loss or damage caused by errors or omissions or by statements resulting from negligence, accident, or any other cause. Brady further assumes no liability arising out of the application or use of any product or system described, herein; nor any liability for incidental or consequential damages arising from the use of this document.

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
WARNUNGEN

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AVISOS

Register your Brady LabPal™ on-line at www.bradylabid.com to receive FREE impact guards and your warranty.

POWER: Six AA 1,5V  Alkaline Batteries or AC Adapter 9V ; 2,2A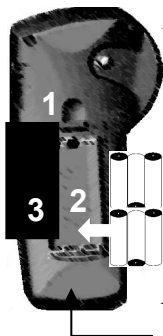

Batteries


1. Open battery compartment.
2. Insert six AA alkaline batteries.
3. Close battery compartment.

-or-

A/C Power


Plug in *optional* A/C Power Supply

DISPLAY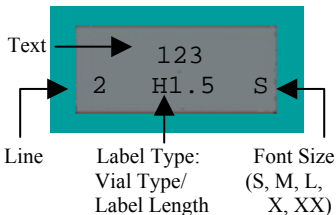

SET TIME & DATE


1. Press  + 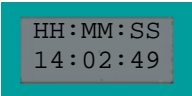

HH : MM : SS
14 : 02 : 49

2. Enter time in military format. e.g.:
5 AM = 05:00:00
1 PM = 13:00:00

3. Press 

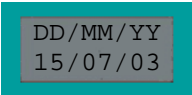DD / MM / YY
15 / 07 / 03

4. Enter date DD/MM/YY

5. Press 

INSERT TIME & DATE

PRESS

 + 

PRINTS

14:02

 + 

15/Jul/03

To lockout access to changing the time/date stamp, press this key combination *simultaneously*: **U, M, Y.**

CREATE LABELS

1. Press

POWER 

2. *Press

LABEL TYPE

Label Types

Vt Gen	Hz Ban	
Vt 1.5ml	Hz 0.6ml	Vial Sizes
Vt 2-4ml	Hz 1.5ml	
	Hz 2-4ml	
Vt 23 mm	Hz 23 mm	Fixed Lengths
Vt 42 mm	Hz 42 mm	
Vt 53 mm	Hz 53 mm	
Vt 66 mm	Hz 66 mm	

3. *Press

FONT SIZE

7 pt (S) 10 pt (M), 14 pt (L),
20 pt (X), 28 pt (XX)

4. Enter Text

Max. Printable Lines = 4

5. Press

Locking Lever must be locked to print. See **Change Label Cartridge.**


6. Cut Label-slide cut lever up

*Press key to scroll through selections.
FUNCTION + key scrolls backwards.

SERIALIZATION


Use arrows ◀ ▶ to place cursor under serial character.

Press **SERIAL** to print each consecutive serialization.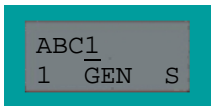

Multiple Serializations


Press **FUNCTION** + **SERIAL**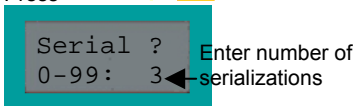

Prints as: ABC1 ABC2 ABC3


Multiple Copies

Press **FUNCTION** + **PRINT**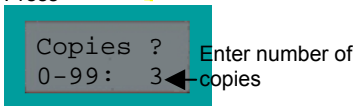

Prints as: ABC4 ABC4 ABC4

CHANGE LABEL CARTRIDGE

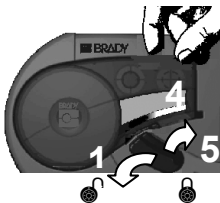1. Unlock Lever
2. Push Release Button (on back)
3. Remove Old Cartridge
4. Insert New Cartridge
5. Lock Lever

KEYPAD

FUNCTION	PRESS
Uppercase	FUNCTION + ▲
Lowercase	FUNCTION + ▼
New Line	ENTER
Space	SPACE
Font Size	FONT SIZE
Delete Character	CLEAR DELETE
Clear Legend	FUNCTION + CLEAR DELETE
Feed	FUNCTION + 1 FEED

SYMBOLS

Press **FUNCTION** + Character

CHAR	SYMB	CHAR	SYMB	CHAR	SYMB
A	%	L	Σ	W	Ö
B	&	M	#	X	Û
C	,	N	/	Y	Æ
D	.	O	*	Z	Ⓢ
E	α	P	(2	²
F	β	Q)	3	↕**
G	γ	R	:	7	-
H	Δ	S	°	8	+
I	λ	T	Ä	9	=
J	μ	U	Å	0	±
K	ρ*	V	Ñ		

*Prints as ρ

** Prints as ↑, ↓, ←, or →. Press key repeatedly to get desired arrow.

GLOBAL CHARACTERS

Use arrows ◀ ▶ to place cursor under character.

Press **FUNCTION** + ⓘ repeatedly until the desired character displays.

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Battery: This product contains a Lithium battery. Dispose of this Lithium battery according to your state or local regulations.

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Caution: Changes or modifications not expressly approved by Brady could void the user's authority to operate the printer.

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US

Note: This printer has been tested and found to comply with the limits for a Class A digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the printer is operated in a commercial environment. This printer generates, uses, and can radiate radio frequency energy and, if not used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this printer in a residential area is likely to cause harmful interference in which case users will be required to correct the interference at their expense.

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Europe

Warning: This is a class A product. In a domestic environment this product may cause radio interference in which case the user may be required to take adequate measures.

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P.O. Box 2131 • Milwaukee, Wisconsin 53201-2131 U.S.A.